


## Bolt Press & Hydraulic Die – *Low-Cost Pellet Preparation*

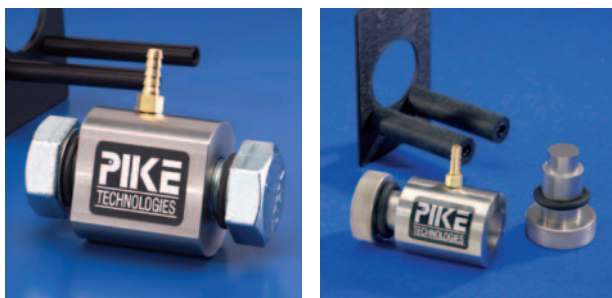

The PIKE Technologies Bolt Press and Hydraulic Die are low-cost tools for making KBr pellets for transmission FTIR analysis.

The press and die consist of a stainless steel barrel with two hardened and polished 13 mm rams. The barrels are equipped with a fitting which allows evacuation of air while the pellet is formed. For the Bolt Press, the pressure is applied to the sample by tightening the bolts against each other with standard 15/16" wrenches – included. For the Hydraulic Die the pressure is applied to the sample by placing in a hydraulic press – up to 10,000 psi. Once a clear pellet is formed, the rams are removed and the sample is analyzed while still in the barrel (barrel is placed directly in the beam using the Press Holder with a standard 2" x 3" slide mount).

Both accessories form a 13 mm pellet. The PIKE Technologies Bolt Press and Hydraulic Die are both equipped with a holder for your FTIR sample compartment.

### ORDERING INFORMATION

#### Pellet Press

PART NUMBER	DESCRIPTION
-------------	-------------

161-2500	Bolt Press for 13 mm pellets
----------	------------------------------

161-3500	Hydraulic Die for 13 mm pellets
----------	---------------------------------

*Notes: The Bolt Press includes evacuable barrel, 2 anvil bolts, 2 15/16" wrenches, and Bolt Press Holder. The Hydraulic Die includes evacuable barrel, 2 rams and Hydraulic Die Holder.*

#### Pellet Press Options and Replacement Parts

PART NUMBER	DESCRIPTION
-------------	-------------

160-8010	KBr Powder, 100 g
----------	-------------------

161-5050	Agate Mortar and Pestle, 50 mm
----------	--------------------------------

161-2511	Replacement 15/16" wrench
----------	---------------------------

161-2520	Bolt Press Holder
----------	-------------------

*Notes: For other options for the Pellet Presses, please contact PIKE Technologies.*

## Hand Press – *for Making Smaller Pellets*



The PIKE Technologies Hand Press is an ideal solution for laboratories that require only occasional preparation of KBr pellets and cannot justify the expense of a hydraulic press.

The Hand Press is an efficient, reliable and inexpensive tool which simplifies making small pellets. It consists of a long stainless steel barrel and movable stage controlled by a lever capable of applying high pressure to the KBr/powder mixture. The Hand Press comes complete with three standard die sets (1, 3 and 7 mm). The pellet preparation involves loading of the powdered sample into the die chamber, placement of the upper anvil in the press and application of hand pressure to the lever (this is sufficient to provide clear, high quality KBr disks). The Die Collar with the formed pellet is removed from the press and in most cases it can be placed directly in the beam of the spectrometer for analysis. The Hand Press is equipped with a platen position dial for adjustment of the force applied to the die and reproducible sample preparation.

### ORDERING INFORMATION

#### Hand Press

PART NUMBER	DESCRIPTION
-------------	-------------

161-1100	Hand Press for 7 mm, 3 mm, and 1 mm pellets <i>Includes: 7 mm, 3 mm, and 1 mm die sets, anvils, die collars, anvil ejectors and Dual Pellet Holder</i>
----------	---

#### Hand Press Options and Replacement Parts

PART NUMBER	DESCRIPTION
-------------	-------------

161-5700	Dual Pellet Holder for 7 mm, 3 mm, and 1 mm pellets
----------	---

161-1018	Single Pellet Holder for 7 mm pellets
----------	---------------------------------------

160-8010	KBr Powder, 100 g
----------	-------------------

161-5050	Agate Mortar and Pestle, 50 mm
----------	--------------------------------

161-1027	Hand Press Body
----------	-----------------

161-1028	1 mm die set
----------	--------------

161-1024	3 mm die set
----------	--------------

161-1010	7 mm die set
----------	--------------

*Notes: For other options for the Hand Press, please contact PIKE Technologies.*